

## PARTICLE SIZE

## ZETA POTENTIAL

MOLECULAR WEIGHT

## RHEOLOGICAL PROPERTIES

# BeNano Series

# Be the Nanoparticle Expert You Need

**Bettersize**  
BETTER PARTICLE SIZE SOLUTIONS

# ADDITIVE

The BeNano Series is widely applied in academic research and manufacturing processes across various fields, including but not limited to pharmaceuticals, nanomaterials, household chemicals, foods and beverages, and academia.



## Pharmaceuticals

**Sample:** Fat emulsions, liposomes, vaccines, hydrogels, etc.

In the field of pharmaceuticals, the BeNano can evaluate systematic stability and alleviate risks in formulations by characterizing size and zeta potential, accelerating the R&D process. The size and size distribution of drugs and drug delivery systems are closely related to the manufacturing process and significantly impact bioavailability, efficacy, and immune response produced by the final product.



## Proteins & Polypeptides

**Sample:** Lysozyme, Human serum albumin, Immunoglobulin G (IgG), etc.


For a protein formulation, it's important to detect subtle changes in size and stability to ensure efficacy and safety. Proteins in poor formulations are likely to form aggregates, which will reduce the efficacy of protein drugs and, even worse, cause immunological reactions and toxicity. To prevent these issues, the BeNano provides rapid access to the characterization of proteins in terms of size and stability information.



## Paints, Inks & Coatings

**Sample:** Oil-based and water-based paints, organic pigments, ceramic inks, etc.

The size and size distribution of paints, inks, and coatings are crucial indicators for ensuring a long shelf life and optimal performance. Poor product quality may lead to aggregation, color inconsistencies, and blockages in channels or nozzles. By using the BeNano to measure these properties, manufacturers can improve formulation development and enhance product performance.

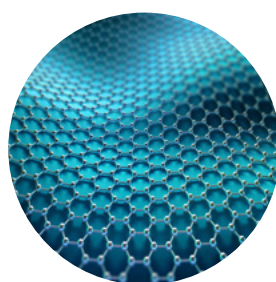

## Nanomaterials

**Sample:** Silver nanoparticles (AgNPs), quantum dots, titanium dioxide, zinc oxide, synthetic silica, etc.

Nanomaterials have wide applications in emerging technologies such as nanoelectronics, nanophotonics, energy conversion, etc. Many of the physicochemical properties associated to nanomaterials are strongly dependent upon the size and particle-particle interaction. With the BeNano, researchers could easily carry out the measurements of size and zeta potential.

# APPLICATIONS



## Household Chemicals

**Sample:** Cosmetics, shampoos, detergents, etc.

Nanomaterials dispersed in sun creams block ultraviolet radiation from the sun. The smaller the particle size, the larger the surface area and the smoother the cream feels. Surfactants in the detergents may remove oil contamination by forming microemulsions. The decontamination effect depends on the type of surfactants used and the size of emulsion droplets. The characterization of nanomaterials is related to all aspects of life.



## Foods and Beverages

**Sample:** Soft drinks, dairy products, confectioneries, plant-based products, etc.

Studies on food and beverages can be performed using the BeNano for characterizing size and zeta potential, in order to optimize the stability of dispersions and emulsifications, which improves the appearance, taste and mouthfeel, and also prolongs the shelf life of products.



## Abrasives

**Sample:** Nano alumina, nano silicon carbide, nano diamond, cubic boron nitride (CBN), etc.

Nano abrasives are extensively used for high-precision polishing and surface finishing of materials such as optical lenses, crystals, gemstones, semiconductors, etc. The stability of polishing slurry is significant for preventing the formation of aggregates that may lead to scratches on workpieces. The BeNano is capable of characterizing the size and zeta potential of slurries, even those with high concentrations.



## Academia

**Sample:** Fundamental and frontier research related to size, zeta potential, molecular weight, etc.

Measurements of size, zeta potential and molecular weight derived from the BeNano provide a powerful tool for academic research, ranging from verifying theory extrapolations to exploring novel synthetic substances. The accurate and highly reproducible data generated by the BeNano ensure the authenticity and reliability of the research results.



# Unlock Greater Research Potential With BeNano

- **Advanced ELS Technology: PALS**

PALS technology can effectively distinguish and extract the electrophoretic behavior even for sample with weak electrophoretic mobilities, either close to isoelectrical point or with high salinity environment.

- **Advanced DLS Technology: Backscattering Detection**

Backscattering DLS optics can detect much larger scattering volume compared to 90-degree optics. Combined with movable measurement position, backscattering DLS offers much higher sensitivity and high turbidity sample measurement capacity.

- **DLS Flow Mode**

Combined with GPC/SEC, or FFF, DLS flow mode can offer a high-resolution size result better than 1.3x to distinguish monomers, dimers and aggregates in a polydisperse sample.

- **Ultra Low Sample Volume Required**

Measuring trace amount of sample is required for early-stage R&D in pharmaceutical industry and academia. With the capillary sizing cell, only 3 to 5  $\mu\text{L}$  of sample is needed for precise size measurement.

- **Temperature Trend Measurement**

For thermal sensitive samples, a temperature trend can be performed easily with a programmed SOP. The BeNano can detect the temperature transition point of the size results, which is the aggregation temperature for protein samples.



Model	Technology		
	90° DLS & SLS	173° DLS & SLS	12° ELS & PALS
BeNano 180 Zeta Pro	✓	✓	✓
BeNano 180 Zeta		✓	✓
BeNano 90 Zeta	✓		✓
BeNano Zeta			✓
BeNano 180 Pro	✓	✓	
BeNano 180		✓	
BeNano 90	✓		

## • Research Level Software

The BeNano software can evaluate and process scattered light signals intelligently to improve the signal quality and result stability. Various built-in calculation modes can cover multiple scientific research and application fields.

## • Stable and Durable Optical Bench

The BeNano adopts a 50mW solid-state laser, a single-mode fiber system and a high-performance APD detector, providing stable, wide-ranging, and highly redundant detection capabilities.

## Measurable Parameters

Hydrodynamic diameter  **$D_H$**

Polydispersity index **PDI**

Intensity, volume, area and number distributions

Diffusion coefficient  **$D$**

Interaction parameter  **$k_D$**

Zeta potential and its distribution

Isoelectric point **IEP**

Molecular weight  **$M_w$**

Viscoelastic modulus  **$G'$ ,  $G''$**



Key Function				
Particle Size	Zeta Potential	Molecular Weight	DLS Microrheology	DLS Flow Mode (Optional)
✓	✓	✓	✓	✓
✓	✓	✓	✓	✓
✓	✓	✓	✓	✓
	✓			
✓		✓	✓	✓
✓		✓	✓	✓
✓		✓	✓	✓



# Particle Size Measured by DLS

Dynamic light scattering (DLS), also referred to as photon correlation spectroscopy (PCS) or quasi-elastic light scattering (QELS), is a technique used to measure Brownian motion in a dispersant. It is based on the principle that smaller particles move faster while larger particles move slower. The scattering intensities of the particles are detected by an avalanche photodiode (APD) and then converted into a correlation function using a correlator. From this correlation function, a mathematic algorithm can be applied to obtain the diffusion coefficient ( $D$ ). The hydrodynamic diameter ( $D_H$ ) and its distribution can be calculated using the Stokes-Einstein equation, which relates the diffusion coefficient to the particle size.


$$D = \frac{K_B T}{3\pi\eta D_H}$$

## Applications


- 1 Particle size and distribution of polymers, colloids, biomacromolecules, etc.
- 2 Research on thermal-sensitive systems, such as PNIPAM polymer.
- 3 Studies on the polymerization process and reaction mechanisms.
- 4 Kinetics analysis of self-assembly, polymerization, and depolymerization, etc.

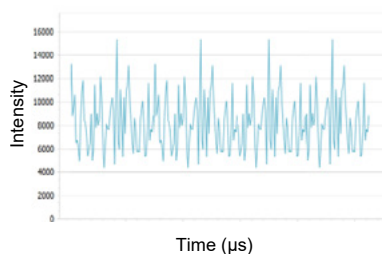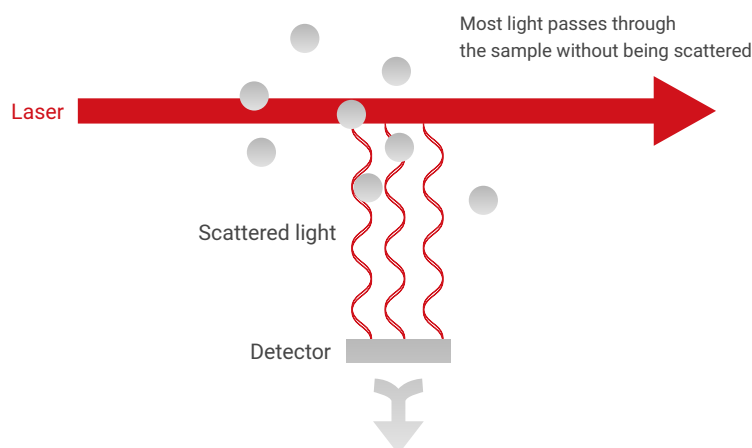


Intensity Fluctuations of Small Particles

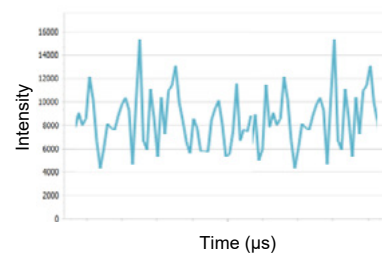

Intensity Fluctuations of Large Particles

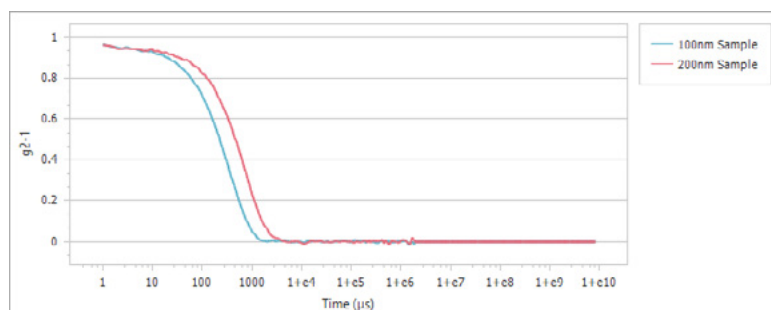

In correlation functions, a 200nm sample decays slower than a 100nm sample due to slower intensity fluctuation.

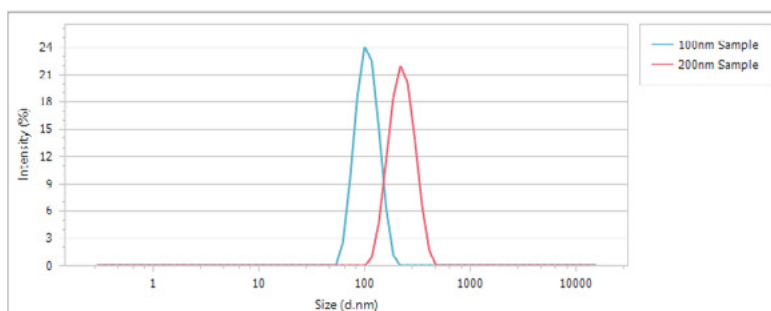

From the above correlation functions, the size and size distribution can be determined using the Stokes-Einstein equation.

# Backscattering Detection Technology

## Features

- **Wider Concentration Range**

By optimizing the detection position, the highly concentrated samples can be detected near the edge of the sample cell, effectively minimizing the multiple light scattering effect.

- **Higher Sensitivity**

8 - 10 times scattering volume and around 10 times sensitivity as compared to the traditional 90° optical design.

- **Higher Size Upper Limit**

It mitigates multiple light scattering from large particles and, to some extent, reduces the number fluctuation of large particles due to the much larger scattering volume.


- **Better Reproducibility**

The DLS backscattering technology is less influenced by dust contaminants and unevenly distributed agglomerates and provides better reproducibility.



## Intelligent Search for the Optimal Detection Position

The software automatically determines the optimal detection position based on the size, concentration, and scattering ability of the sample to achieve the highest measurement accuracy and offer flexibility in detecting different types and concentrations of samples. This feature is particularly useful when dealing with a variety of samples, each with its unique scattering properties and concentrations.

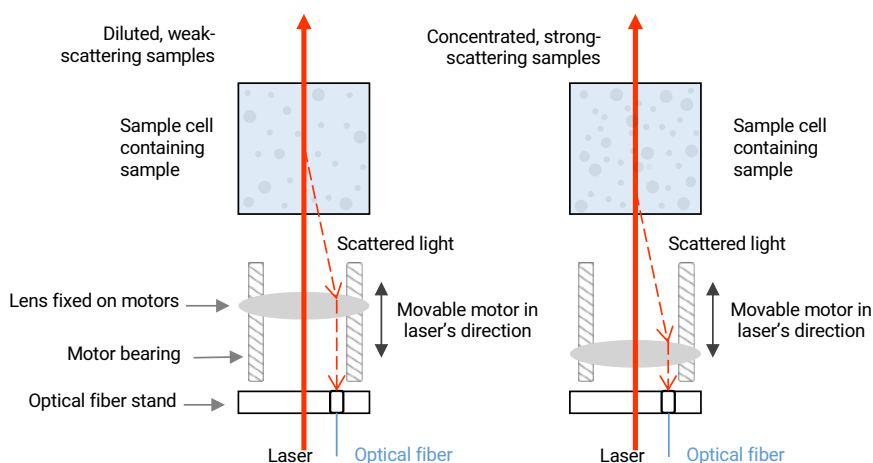

- 1 **The detection point in the middle of the sample cell**

This leads to a large scattering volume that increases instrument sensitivity and is suitable for detecting dilute samples with weaker scattering effects.

- 2 **The detection point at the edge of the sample cell**

This avoids the multiple scattering effect of high concentration samples, ensuring accurate and repeatable particle size results.


# DLS Flow Mode

DLS flow mode provides a high-resolution size result of a complex, polydisperse system. When combined with front-end separation equipment such as GPC/SEC or FFF, particles are separated into monodisperse fractions and flow through the BeNano in sequence by size. The size of each fraction is continuously measured and summed into a high-resolution size distribution.

BeNano can acquire RI or UV signals, offering a more accurate volume and number distributions independent of algorithm compared to a batch-mode measurement.




Integrated use of BeNano and SEC for particle characterization

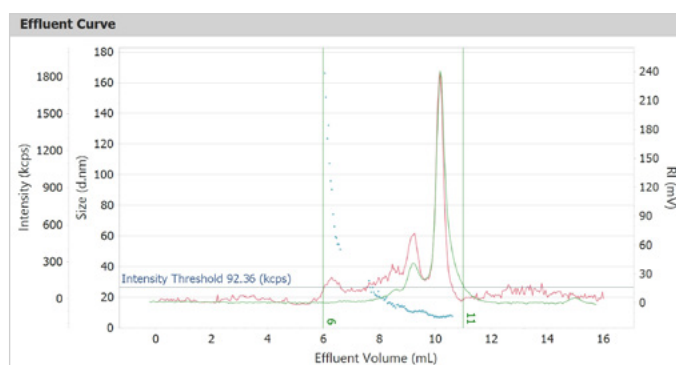

Effluent curves of size, intensity, and refractive index (RI)

## Applications

- 1 Characterizing particulates and polymers on size and dispersity
- 2 Distinguishing monomers, dimers and aggregates of proteins

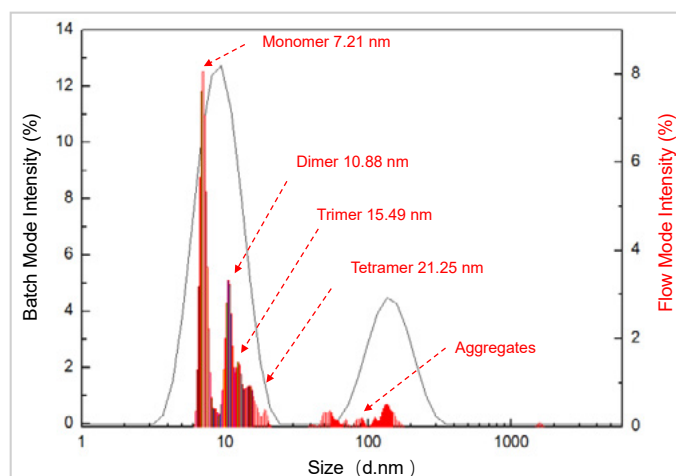


Flow mode provides higher resolution for characterizing BSA on size and distribution compared with batch mode




# High-resolution Size Measurement

## Features


- DLS analyzer connecting with GPC/SEC, FFF, etc.
- Receiving up to 3 signals from RI, UV or other detectors
- 27uL low-volume flow cell to avoid band broadening
- Size resolution as high as 1.3 : 1
- Size distributions weighted by number and volume besides intensity
- Suitable for complex, polydisperse systems such as proteins, polymers, etc.

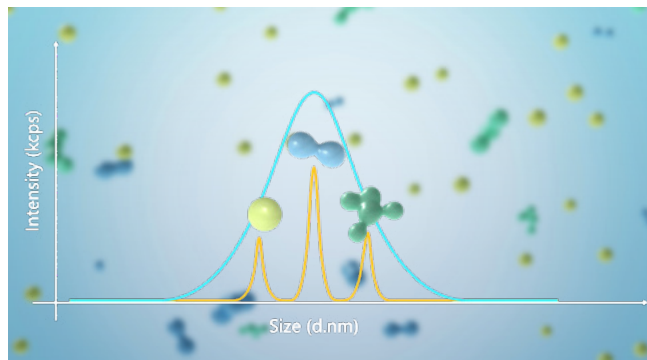

High-resolution size distribution achieved through flow mode

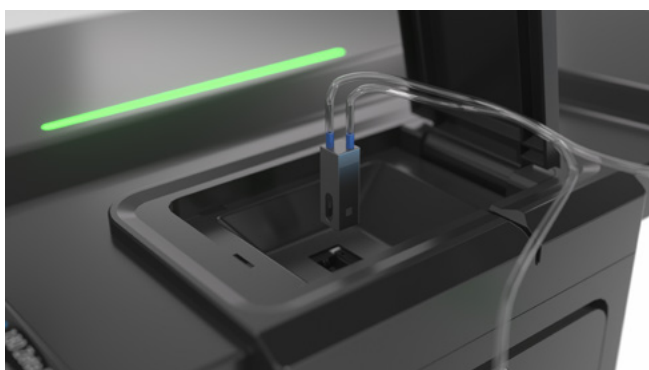

Particles remain in flow during measurement

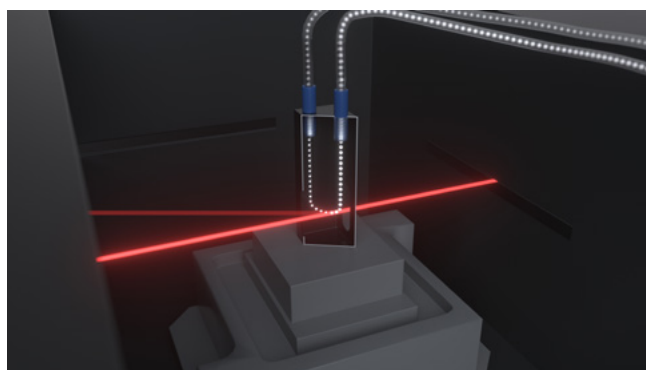


Following separation, particle sizes are determined

## Report Designer

The built-in report designer provides a range of editable templates, enabling users to customize reports to their needs.


- 30 report templates cover all measurement results
- Flexible editing, one-click saving, and easy creation of user-defined reports
- Customizable report logo

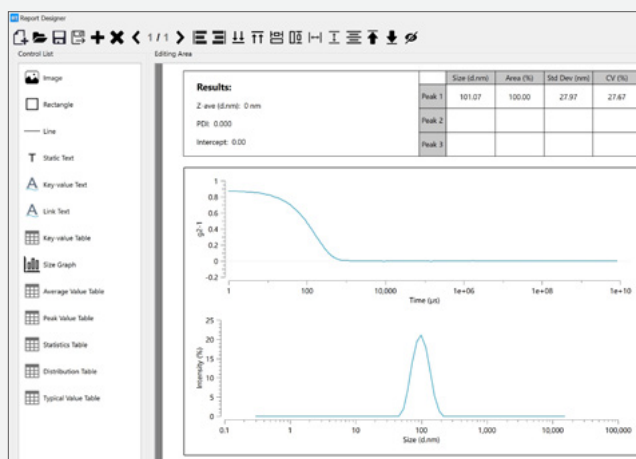


# Zeta Potential Measured by ELS

In aqueous systems, charged particles are surrounded by counterions that form an inner Stern layer and an outer shear layer. Zeta potential is the electrical potential at the interface of the shear layer. A higher zeta potential indicates greater stability and less aggregation of the suspension system. Electrophoretic light scattering (ELS) measures electrophoretic mobility via Doppler shifts of scattered light, which can be used to determine the zeta potential of a sample by Henry's equation.

$$\mu = \frac{2\varepsilon_r\varepsilon_0\zeta}{3\eta} f(k\alpha)$$

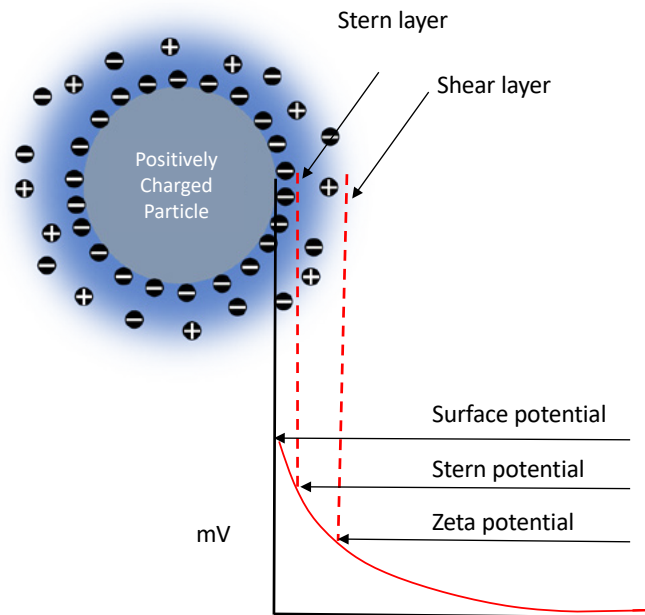

Potential Distribution at Particle Surface

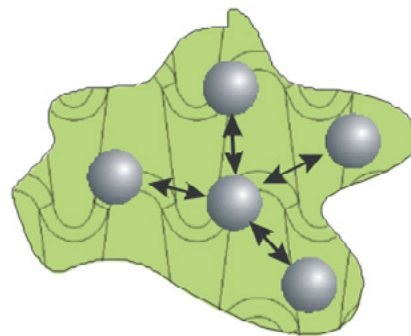


Electrostatic Forces Between Particles

## Applications

- 1 Particle suspensions, including colloids, proteins, nanometals, etc
- 2 Industries such as chemicals, biology, water treatment, paints, etc
- 3 Product stability control and monitoring
- 4 Stability research and control of suspension system
- 5 Surface property and modification study

## Colloidal Stability

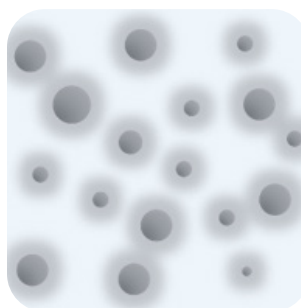

**Stable particle system**

- High repulsion force of particles
- High zeta potential



**Unstable particle system**



- Flocculation, aggregation, sedimentation
- Low or zero zeta potential

# PALS: Phase Analysis Light Scattering

Phase analysis light scattering (PALS) is an advanced technology based on the traditional ELS technology, which has been further developed by Bettersize to measure the zeta potential and its distribution of a sample.

## Features & Benefits

- Accurate measurement of samples with low electrophoretic mobility
- Effective for samples in organic solvents with low dielectric constant
- More accurate results for samples with high conductivity
- Effectively measures the zeta potential of particles whose charge approaches the isoelectric point

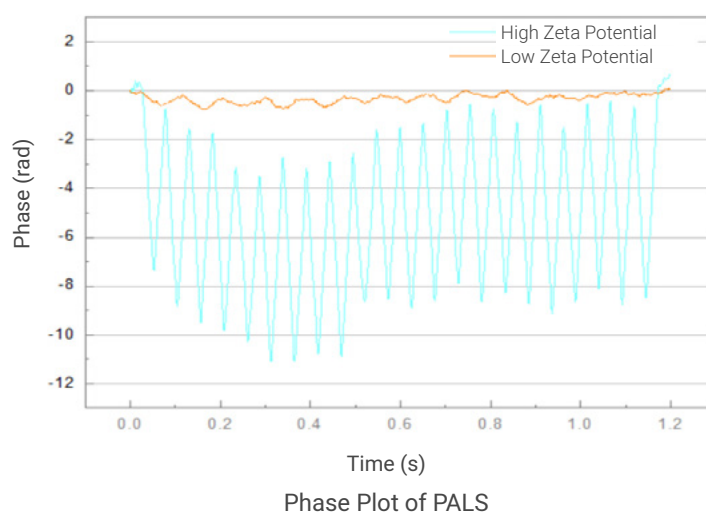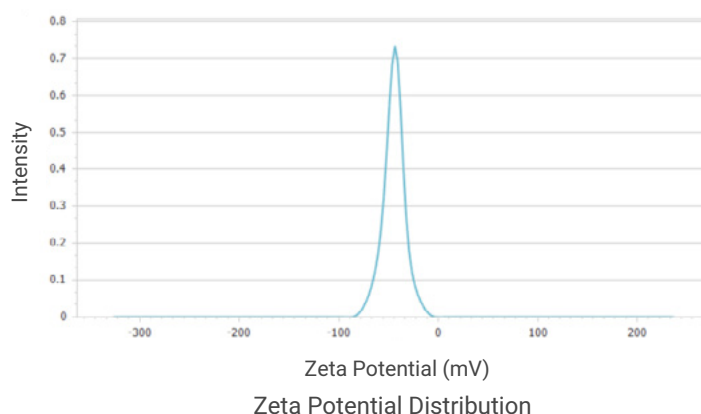


# Static Light Scattering (SLS)

Static light scattering (SLS) is a technology that measures the scattering intensities, weight-average molecular weight ( $M_w$ ) and second virial coefficient ( $A_2$ ) of the sample through Rayleigh equation:

$$\frac{Kc}{R_\theta} = \frac{1}{M_w} + 2A_2c$$


where  $c$  is the sample concentration,  $\theta$  is the detection angle,  $R_\theta$  is the Rayleigh ratio used to characterize the intensity ratio between the scattered light and the incident light at the angle of  $\theta$ ,  $M_w$  is the sample's weight-average molecular weight,  $A_2$  is the second virial coefficient, and  $K$  is a constant related to  $(dn/dc)^2$ .

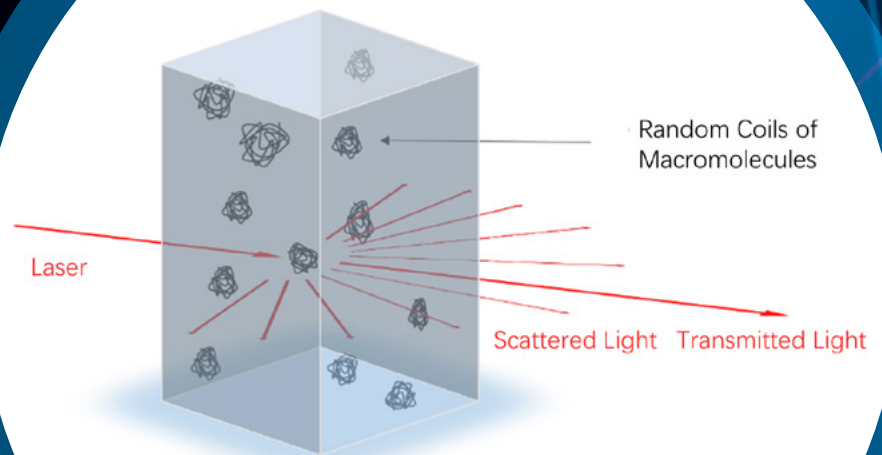

Scattered Light of Macromolecules

## Applications

- 1 **Chemical engineering:**  
characterization of polymers, micelles and macromolecules
- 2 **Life science:**  
characterization of proteins, polypeptides, and polysaccharides
- 3 **Pharmaceuticals:**  
research on aggregation and stability of drugs and biomacromolecules

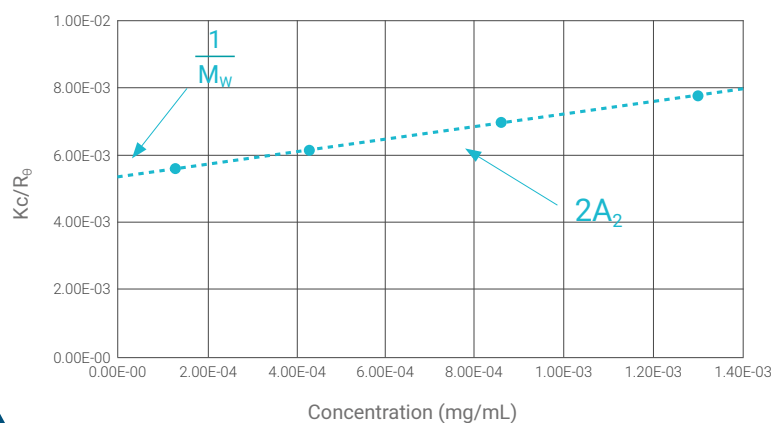

Debye Plot

# Microrheology Measured by DLS

## Dynamic Light Scattering

Microrheology (DLS Microrheology) is an economical and efficient technique that utilizes dynamic light scattering to determine rheological properties. By analyzing the Brownian motion of colloidal tracer particles, information about the viscoelastic properties of the system, such as viscoelastic modulus, complex viscosity and creep compliance, can be obtained with the generalized Stokes-Einstein equation.

## Features & Benefits

- Investigates rheological behaviors by measuring the thermally-driven motion of tracer particles within a material being studied
- Facilitates the measurement across a wide range of frequencies
- Applies low stress to tracer particles
- Requires only a microliter-scale sample volume
- Complements mechanical rheology results
- Suitable for weakly-structured samples

## Applications

- 1 Polymer solution
- 2 Protein solution
- 3 Gel system

## Stokes-Einstein equation

$$G^*(\omega) = \frac{k_B T}{\pi R i \langle \Delta r^2(i\omega) \rangle} = G'(\omega) + iG''(\omega)$$


## Microrheology Information

Elastic (storage) modulus  $G'$


Complex viscosity  $\eta^*$

Viscous (loss) modulus  $G''$


Creep compliance  $J$

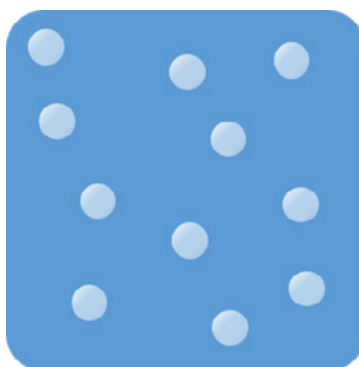


Particles in pure viscous fluids

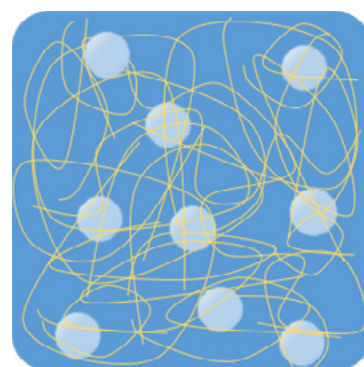

Particles in complex viscoelastic fluids

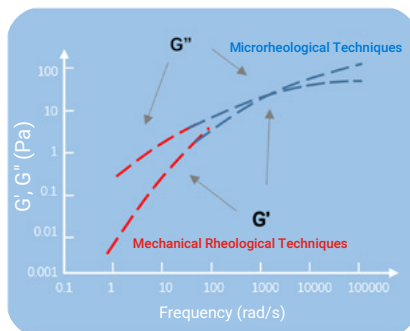

$G'$  and  $G''$  curves of complex fluids

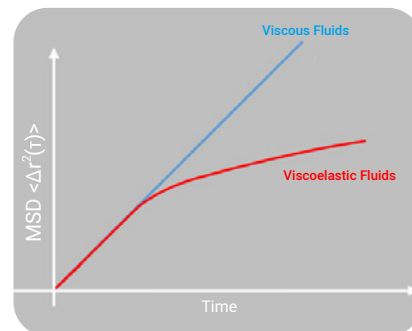

MSD curves of particles in pure viscous fluids and complex viscoelastic fluids




# Temperature Trend Measurement

## Measurement Parameters

- Size vs. Temperature
- Zeta Potential vs. Temperature


## Features

- Temperature trend measurement function (-15°C to 110°C)
- Significant for investigating particle size and zeta potential under different temperatures

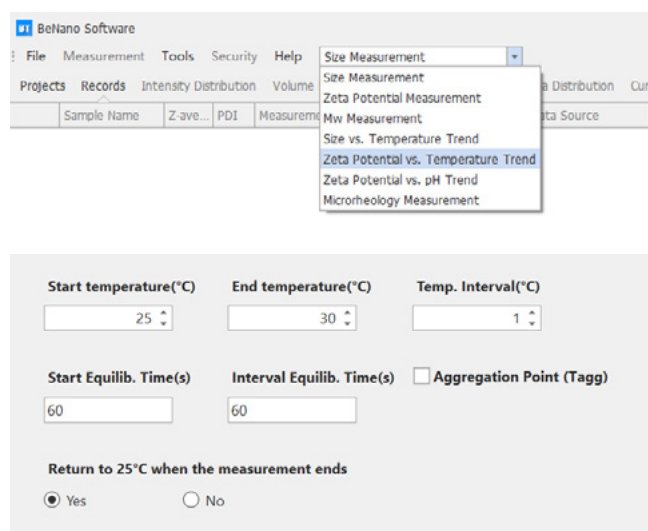

## Benefits


- Easy examination of protein formulation stability
- Accelerates real-time aging through elevated temperature simulation

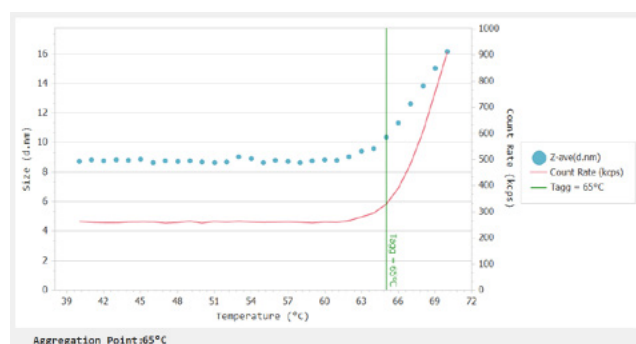

Size vs. Temperature trend measurement of the BSA protein

# A Research Level Software

- SOP guarantees measurement accuracy and completeness
- Automatic calculation of mean and standard deviation for results and statistics
- Comparison of results from multiple runs through statistics and overlay functions
- Real-time display of information and results
- Over 100 available parameters that meet research, QA, QC, and production needs
- Free lifelong upgrades provided

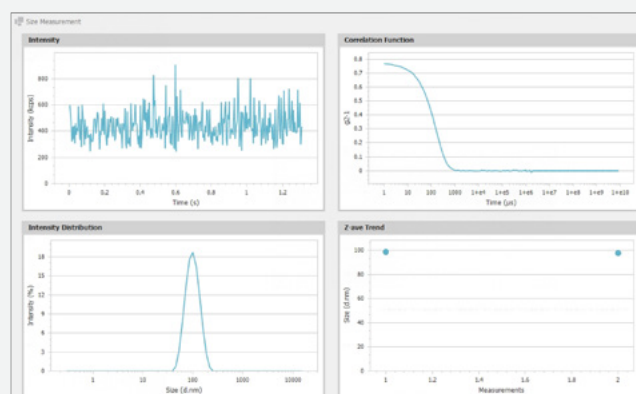

# pH Trend Measurement

## Measurement Parameters

- Zeta potential vs. pH
- Isoelectric point
- Conductivity vs. pH

## Features

- High-precision ternary titration pumps
- Controllable peristaltic pump with high flow capacity and high flow rate
- General-purpose electrode
- Automated titrant selection based on initial and target pH using intelligent software


Start pH: 3, End pH: 2, ☒ Titrate from current pH

☒ pH Interval: 1, ☐ pH Points: 5

pH Tolerance: 0.2

## Benefits

- Completes measurements within a shorter time
- Improves consistency and repeatability of results
- Reduces the workload of researchers
- Simplifies qualifications needed for operators
- Reduces exposure to corrosive liquids

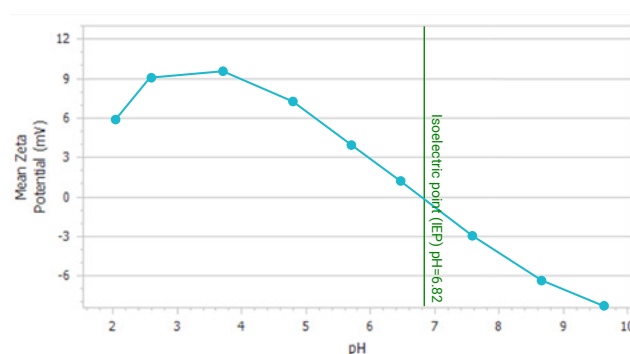

## Compliance With FDA 21 CFR Part 11

The BeNano software system is compliant with 21 CFR Part 11 regulations, which restricts access to authorized individuals through a username and password system for electronic record signing, access logs, change logs, or operation execution. An activation code can be used to upgrade security settings and ensure compliance, and an "audit trail" can be viewed to ensure proper management and maintenance of system security and data integrity.

Activate 21 CFR feature

Status: Security Activated

Instrument Serial No.  
02 . [ ] . [ ] . [ ]

License Key  
[ ]

Activate Security



# CELLS

## Standard



### Folded Capillary Cell BT-C1

For aqueous samples in zeta potential measurement

Material: PC, Gold-plated Phosphor Bronze

Sample Volume: 0.75 mL

Temperature Range: -15 - 70 °C



### PS Cuvette

For aqueous samples in size measurement

Material: Polystyrene

Sample Volume: 1 - 1.5 mL

Temperature Range: -15 - 70 °C



### Glass Cuvette

For aqueous and organic samples in size measurement

Material: Quartz

Sample Volume: 1 - 1.5 mL

Temperature Range: -15 - 110 °C



### Capillary Sizing Cell


For aqueous and organic samples with ultra-micro volume required in size measurement

Material: Quartz

Sample Volume: 3 - 5  $\mu$ L

Temperature Range: -15 - 70 °C

## Optional

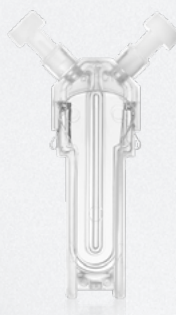

### Folded Capillary Cell BT-C1-Pt

For aqueous samples in zeta potential, high salinity tolerance

Material: PC, Platinum

Sample Volume: 0.75 mL

Temperature Range: -15 - 70 °C




### Micro-volume Cuvette

For aqueous samples with micro volume required in size measurement

Material: PMMA

Sample Volume: 40 - 50  $\mu$ L

Temperature Range: -15 - 70 °C

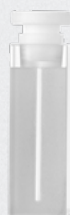

### Micro-volume Glass Cuvette

For aqueous and organic samples with micro volume required in size measurement

Material: Quartz

Sample Volume: 25  $\mu$ L

Temperature Range: -15 - 110 °C



### Dip Cell

For aqueous and organic samples in zeta potential measurement

Material: PEEK, Platinum

Sample Volume: 1 - 1.5 mL

Temperature Range: -15 - 70 °C



# BAT-1 AUTOTITRATOR

The BAT-1 autotitrator, a highly recommended accessory for the BeNano series, is an ideal option for analyzing the zeta potential as a function of pH and determining the isoelectric point (IEP) of a colloidal system. It is equipped with three high-precision titration pumps with a precision of 0.28  $\mu\text{L}$ , and the system can select the optimal titrant and adjust the addition volume intelligently to optimize titration efficiency and accuracy.

**Making pH titration more efficient, more accurate and safer** 



## Flow Cell

For aqueous and organic samples in DLS flow mode measurement

Material: Quartz

Sample Volume: 27  $\mu\text{L}$

Temperature Range: 0 - 70  $^{\circ}\text{C}$



## Micro-volume Fluorescence-filtered Cell

For aqueous samples in size measurement

Material: Quartz

Sample Volume: 16  $\mu\text{L}$

Temperature Range: 0 - 90  $^{\circ}\text{C}$

Measurement Angle: 90 $^{\circ}$ , 173 $^{\circ}$



## Micro-volume VH Cell

For aqueous samples in size measurement

Eliminating the vertical element of scattered light

Material: Quartz

Sample Volume: 16  $\mu\text{L}$

Temperature Range: 0 - 90  $^{\circ}\text{C}$

Measurement Angle: 90 $^{\circ}$



## Micro-volume VV Cell

For aqueous samples in size measurement

Eliminating the horizontal element of scattered light

Material: Quartz

Sample Volume: 16  $\mu\text{L}$

Temperature Range: 0 - 90  $^{\circ}\text{C}$

Measurement Angle: 90 $^{\circ}$ , 173 $^{\circ}$

Prepare the titrants



Calibrate the pH probe



Create an SOP



Start the test and wait



Check the results



Function	Parameter	BeNano 180 Zeta Pro	BeNano 180 Zeta	BeNano 90 Zeta
Size measurement	Size measurement range	0.3 nm - 15 µm*	0.3 nm - 10 µm*	0.3 nm - 15 µm*
	Sample volume	3 µL - 1 mL*	40 µL - 1 mL*	3 µL - 1 mL*
	Detection angle	90° & 173° & 12°	173° & 12°	90° & 12°
	Analysis algorithm	Cumulants, General Mode, CONTIN, NNLS	Cumulants, General Mode, CONTIN, NNLS	Cumulants, General Mode, CONTIN, NNLS
	Upper limit of concentration range	40% w/v*	40% w/v*	Optically clear†
	Detection position	Movable	Movable	Fixed
Zeta potential measurement	Detection angle	12°	12°	12°
	Zeta potential measurement range	No actual limitation	No actual limitation	No actual limitation
	Electrophoretic mobility	> ± 20 µm·cm/V·s	> ± 20 µm·cm/V·s	> ± 20 µm·cm/V·s
	Conductivity	0 - 260 mS/cm	0 - 260 mS/cm	0 - 260 mS/cm
	Sample volume	0.75 - 1 mL	0.75 - 1 mL	0.75 - 1 mL
	Sample size	2 nm - 120 µm*	2 nm - 120 µm*	2 nm - 120 µm*
Other measurements	Molecular weight (Mw)	342 Da - 2 x 10 <sup>7</sup> Da*	342 Da - 2 x 10 <sup>7</sup> Da*	342 Da - 2 x 10 <sup>7</sup> Da*
	DLS Microrheology	MSD, G', G'', η*, J	MSD, G', G'', η*, J	MSD, G', G'', η*, J
	Viscosity	0.01 cp - 100 cp*	0.01 cp - 100 cp*	0.01 cp - 100 cp*
	Interaction parameter k <sub>D</sub>	No actual limitation	No actual limitation	No actual limitation
	Trend measurement	Time, temperature and pH	Time, temperature and pH	Time, temperature and pH
System parameter	Temperature control range	-15°C - 110°C , ± 0.1°C	-15°C - 110°C , ± 0.1°C	-15°C - 110°C , ± 0.1°C
	Condensation control	Dry air or nitrogen	Dry air or nitrogen	Dry air or nitrogen
	Laser source	50 mW Solid-state laser, 671 nm*, Class 1	50 mW Solid-state laser, 671 nm*, Class 1	50 mW Solid-state laser, 671 nm*, Class 1
	Correlator	Up to 4000 channels, 10 <sup>11</sup> linear dynamic	Up to 4000 channels, 10 <sup>11</sup> linear dynamic range	Up to 4000 channels, 10 <sup>11</sup> linear dynamic range
	Detector	Avalanche photodiode (APD)	Avalanche photodiode (APD)	Avalanche photodiode (APD)
	Intensity control	0.0001% - 100%, manual or automatic	0.0001% - 100%, manual or automatic	0.0001% - 100%, manual or automatic
	Dimensions (L x W x H)	62.5 x 40 x 24.5 cm (26 kg)	62.5 x 40 x 24.5 cm (26 kg)	62.5 x 40 x 24.5 cm (26 kg)
	Power supply	AC 100 - 240 V, 50 - 60 Hz, 4A	AC 100 - 240 V, 50 - 60 Hz, 4A	AC 100 - 240 V, 50 - 60 Hz, 4A
	Compliance	21 CFR Part 11, ISO 13321, ISO 22412, ISO 13099	21 CFR Part 11, ISO 13321, ISO 22412, ISO 13099	21 CFR Part 11, ISO 13321, ISO 22412, ISO 13099

\* Dependent on samples and accessories

† Up to 40% w/v using capillary sizing cell

\* 10mW 633 nm HeNe laser available on request



BeNano Zeta	BeNano 180 Pro	BeNano 180	BeNano 90
N/A	0.3 nm - 15 µm*	0.3 nm - 10 µm*	0.3 nm - 15 µm*
N/A	3 µL - 1 mL*	40 µL - 1 mL*	3 µL - 1 mL*
N/A	90° & 173°	173°	90°
N/A	Cumulants, General Mode, CONTIN, NNLS	Cumulants, General Mode, CONTIN, NNLS	Cumulants, General Mode, CONTIN, NNLS
N/A	40% w/v*	40% w/v*	Optically clear†
N/A	Movable	Movable	Fixed
12°	N/A	N/A	N/A
No actual limitation	N/A	N/A	N/A
> ± 20 µm·cm/V·s	N/A	N/A	N/A
0 - 260 mS/cm	N/A	N/A	N/A
0.75 - 1 mL	N/A	N/A	N/A
2 nm - 120 µm*	N/A	N/A	N/A
N/A	342 Da - 2 x 10 <sup>7</sup> Da*	342 Da - 2 x 10 <sup>7</sup> Da*	342 Da - 2 x 10 <sup>7</sup> Da*
N/A	MSD, G', G'', η*, J	MSD, G', G'', η*, J	MSD, G', G'', η*, J
N/A	0.01 cp - 100 cp*	0.01 cp - 100 cp*	0.01 cp - 100 cp*
N/A	No actual limitation	No actual limitation	No actual limitation
Time, temperature and pH	Time, temperature and pH	Time, temperature and pH	Time, temperature and pH
-15°C - 110°C , ± 0.1°C	-15°C - 110°C , ± 0.1°C	-15°C - 110°C , ± 0.1°C	-15°C - 110°C , ± 0.1°C
Dry air or nitrogen	Dry air or nitrogen	Dry air or nitrogen	Dry air or nitrogen
50 mW Solid-state laser, 671 nm*, Class 1	50 mW Solid-state laser, 671 nm*, Class 1	50 mW Solid-state laser, 671 nm*, Class 1	50 mW Solid-state laser, 671 nm*, Class 1
Up to 4000 channels, 10 <sup>11</sup> linear dynamic range	Up to 4000 channels, 10 <sup>11</sup> linear dynamic range	Up to 4000 channels, 10 <sup>11</sup> linear dynamic range	Up to 4000 channels, 10 <sup>11</sup> linear dynamic range
Avalanche photodiode (APD)	Avalanche photodiode (APD)	Avalanche photodiode (APD)	Avalanche photodiode (APD)
0.0001% - 100%, manual or automatic	0.0001% - 100%, manual or automatic	0.0001% - 100%, manual or automatic	0.0001% - 100%, manual or automatic
62.5 x 40 x 24.5 cm (26 kg)	62.5 x 40 x 24.5 cm (26 kg)	62.5 x 40 x 24.5 cm (26 kg)	62.5 x 40 x 24.5 cm (26 kg)
AC 100 - 240 V, 50 - 60 Hz, 4A	AC 100 - 240 V, 50 - 60 Hz, 4A	AC 100 - 240 V, 50 - 60 Hz, 4A	AC 100 - 240 V, 50 - 60 Hz, 4A
21 CFR Part 11, ISO 13099	21 CFR Part 11, ISO 13321, ISO 22412	21 CFR Part 11, ISO 13321, ISO 22412	21 CFR Part 11, ISO 13321, ISO 22412



[www.bettersizeinstruments.com](http://www.bettersizeinstruments.com)  
[info@bettersize.com](mailto:info@bettersize.com)

**Bettersize Instruments Ltd.**

**Address:** No. 9, Ganquan Road, Jinquan Industrial Park,

Dandong, Liaoning, China

**Postcode:** 118009

**Tel:** +86-755-26926582

**Fax:** +86-415-6170645

**Bettersize Inc.**

**Address:** Suite K-2, 3188 Airway Ave, Costa Mesa, CA 92626,

United States

**Tel:** +1 833-699-7493 (SIZE)

Visit Our BeNano Series Site:



Visit Our Official Youtube Channel:



**Disclaimer:** By using or accessing the brochure, you agree with the Disclaimer without any qualification or limitation. Diligent care has been used to ensure that the information in this brochure is accurate, Bettersize Instruments Ltd. shall not be liable for errors contained herein or for damages in connection with the use of this material. The information on this brochure is presented as general information and no representation or warranty is expressly or impliedly given as to its accuracy, completeness or correctness. It does not constitute part of a legal offer or contract. Bettersize Instruments Ltd. reserves the right to modify, alter, add and delete the content outlined in the brochure without prior notice and without any subsequent liability to the company.

Copyright: © 2024 Bettersize Instruments Ltd. | All Rights Reserved

RIFERIMENTO PER L'ITALIA



**Qi srl**  
t +39 06 9105461  
[www.qitech.it](http://www.qitech.it) | [SalesQi@qitech.it](mailto:SalesQi@qitech.it)